

Quarton inc.

QLM-1125 / QLM-1165 / QLM-1225 User Manual



QLM-1125 / QLM-1165 / QLM-1225

Laser Mounts are compatible with laser module / laser pointer / flashlight with OD 6mm to 22.5mm, it's a ideal choose for mounted onto optical tables.

| Model name | QLM-1125 | QLM-1165 | QLM-1225 |
|------------|--------------|--------------|--------------|
| Dimension | 69*66.4 mm | 69*67 mm | 90*75 mm |
| Weight | 151.4 g | 151.6 g | 168 g |
| OD | Up to 12.5mm | Up to 16.5mm | Up to 22.5mm |




FEATURES

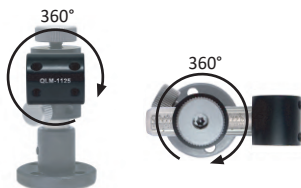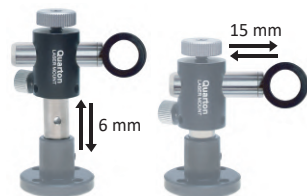

4 directions adjustable.



Can be mounted onto the optical table with screws.
(Screws not included)

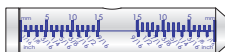


Engraved scale for precision adjustable.

CONTENTS




| NO. | Part Name | QTY | Part Stock Number |
|-----|--------------------------|-----|---|
| 1 | Mount holder | 1 | 405-10008G - 1125 405-10007G - 1165 405-10009G - 1225 |
| 2 | Plastic Screw | 4 | 300246901G |
| 3 | Mount main body | 1 | 405-10006G - 1125&1165 405-10010G - 1225 |
| 4 | Wrench (M2) | 1 | 428-020048-011G |
| 5 | Hard Cover Carrying Case | 1 | 492-5B103G |
| 6 | User Manual | 1 | 494-10709G |

MOUNTING

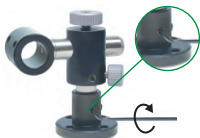

1. Move to desired height and rotate to desired direction, then use wrench tight screws to fix pole firm.

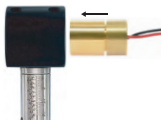

2. Place your laser module/ laser pointer/flashlight to suitable position.



3. Tight screws to fix laser module/laser pointer/ flashlight sturdy.



4. Loose the fixing ring. Rotate the adjustable laser holder to find the angle you want.



5. Tighten the fixing ring to fix the laser holder.

APPLICATION



Flashlight with QLM-1165



Laser pointer with QLM-1165



Ø12.5 with QLM-1165




Ø16 with QLM-1165



Ø16 with QLM-1165



Ø11 with QLM-1125

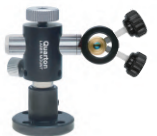

Ø10.5 with QLM-1125



Ø6.5 with QLM-1125