

# Quarton inc.

## QLM-1125 / QLM-1165 / QLM-1225 User Manual


### QLM-1125 / QLM-1165 / QLM-1225

Laser Mounts are compatible with laser module / laser pointer / flashlight with OD 4mm to 22.5mm, it's a ideal choose for mounted onto optical tables.


Model name	QLM-1125	QLM-1165	QLM-1225
Dimension	69*66.4 mm	69*67 mm	90*75 mm
Weight	151.4 g	151.6 g	168 g
OD	Up to 12.5mm	Up to 16.5mm	Up to 22.5mm




### FEATURES

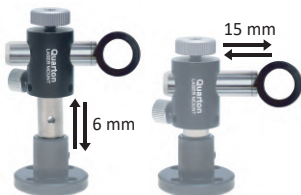

Can be mounted onto the optical table with screws.  
(Screws not included)

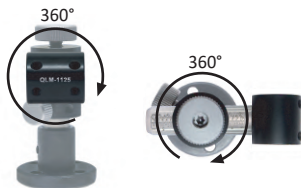

4 directions adjustable.

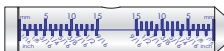


Engraved scale for precision adjustable.

### CONTENTS




NO.	Part Name	QTY	Part Stock Number
1	Mount holder	1	405-10008G
2	Plastic Screw	4	424-40001G
3	Mount main body	1	405-10006G
4	Wrench (M2)	1	428-020048-011G
5	Hard Cover Carrying Case	1	492-5B103G
6	User Manual	1	494-10709G


## MOUNTING

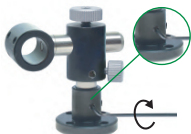


1. Move to desired height and rotate to desired direction, then use wrench tight screws to fix pole firm.

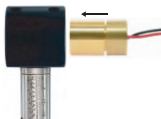

2. Place your laser module/ laser pointer/flashlight to suitable position.

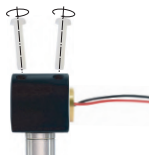

3. Tight screws to fix laser module/laser pointer/ flashlight sturdy.

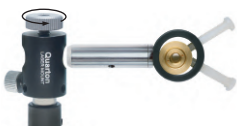

4. Loose the fixing ring. Rotate the adjustable laser holder to find the angle you want.



5. Tighten the fixing ring to fix the laser holder.

## APPLICATION



Flashlight with QLM-1165



Laser pointer with QLM-1165




Ø12.5 with QLM-1165



Ø16 with QLM-1165



Ø16 with QLM-1165

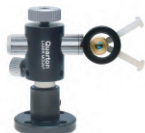

Ø10.5 with QLM-1125



Ø6.5 with QLM-1125



Ø4 with QLM-1125